

Guidelines for collecting and sending tissues (HU,MK)

Guidelines for collecting and sending tissues

All procedures that require aseptic technique

1. Complete soak skin with 70% EtOH;
2. Remove tissues (3-10 gram of skin, pancreas, fat, muscle, Lung, Heart, etc; 5-10 cm vascular tissues of artery, vein, aorta etc. as requested) with a sterile scissors and tweezers.
3. Place tissues in a 50-100 ml sterile tube/bottle containing 40-100 ml of cold DMEM with Gentamycin 200 ug/ml.
4. Seal a tube with parafilm. Gently agitate. Keep a tube containing tissues in cold medium on ICE.
5. Have the separated bottle/tubes (100-200 ml) for different tissues, such as livers, skins or bones.
6. We can provide (1) medium with antibiotics, (2) tubes, and (3) 70% EtOH for use of tissue collection. Please let us know the shipping address.
7. Please note that the organs or tissues CANNOT be frozen during shipment.
8. AT THE SAME DAY send tissues by FedEx Priority Overnight mail.

Shipping Address

Cell Biologics, Inc.
2201 West Campbell Park Drive,
Suite 326
Chicago IL 60612 USA
Phone: 312-226-8198 or 312-981-9263